

SamplerSight Pharma

Batch Sampling Software

Easy to use software for data collection, analysis and reporting

Without measurement there is no control®

SamplerSight Pharma is validatable, user-friendly, Windows-based software that controls syringe operated samplers and collects data from a LiQuilaz® II particle counter. SamplerSight Pharma allows operators to manage sampling requirements for batch-based operations. It also provides a comprehensive view of the batch information with histogram, time plot and tabular data presented in an easy to use format.

Database filtering functions are based on batch, lot and location, allowing a fast response to database queries and sample reporting. Multiple pre-configured or user-defined sample configurations reduce set-up time between batches of different types.

USP 788 sensor calibration is easily accomplished in 20 to 30 minutes. SamplerSight Pharma Software is fully 21 CFR Part 11 compliant, and includes IQ/OQ/PQ documentation. Optional installation setup, validation and training available.



BENEFITS

Cost Reduction

- Dedicated application for batch monitoring reduces integration time and speeds up reporting
- Multi-platform compatibility allows migration to new computer systems, improving reliability

Easy to Use

- Simple-to-use icons make software operation trouble-free
- Reporting tools are predefined to speed up batch reports; user configurable formats allow modifications to fit site demands
- Database filtering functions for batch/lot/location allow fast response to database queries and sample reporting

Increased Productivity

- Real-time data displays allow for a rapid response to particle problems
- Multiple user-defined sample configurations reduce set up time between batches of different types
- A single application supports a variety of products and reduces training time for operators

FEATURES

- Multiple data displays: histogram, time plot, and tabular
- Sample-specific setup configuration, recipes and storage
- User-configurable identification labels
- Batch and filters for data retrieval and reporting
- Comprehensive reporting function with CSV and PDF file export
- Auto-print for each sample
- Manual control diagnostics screen
- Alarm specification and generation functionality

APPLICATIONS

Collecting, storing, and viewing data for:


- Injectable product release
- Liquid production batch monitoring
- Particulate matter in ophthalmic solutions monitoring
- Medical device cleanliness monitoring

SamplerSight Pharma

Batch Sampling Software

Specifications

Sampling systems supported	APSS-2000 with LiQuilaz® II liquid particle counters
On-line RS-485 sensors supported	LiQuilaz II (E-Series and S-Series)
Minimum computer requirements	Please refer to the Microsoft® operating system minimum requirements. RS-232 serial communication port (when used in conjunction with non-Ethernet devices), network interface port
Software required	Windows® 10 or Windows 11 with Microsoft SQL Server 2016 installed (SQL Server Express included with the software)

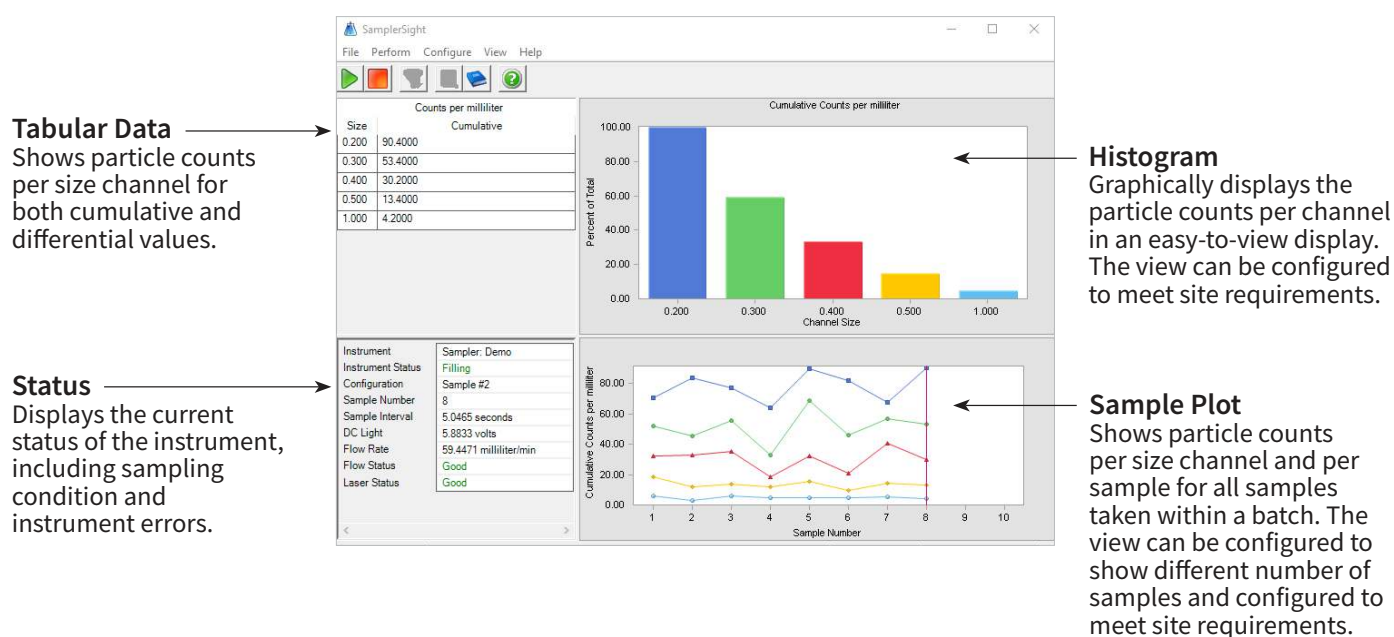

HEADQUARTERS

Boulder, Colorado 80301 USA
T: +1 303 443 7100

Instrument Service & Support
T: +1 303-443-7100 ext 447

Service and Calibration
E: workorders@pmeasuring.com

Technical Support
E: techsupport@pmeasuring.com

Software or System Support
E: ssupport@pmeasuring.com

Customer Service and Order Processing
E: customerorders@pmeasuring.com

www.pmeasuring.com
info@pmeasuring.com

GLOBAL OFFICES

AUSTRIA
T: +43 171 728 285
E: pmsaustria@pmeasuring.com

BELGIUM/BENELUX
T: +32 10 23 71 56
E: pmsbelgium@pmeasuring.com

BRAZIL
T: +55 11 5188 8227
E: pmsbrazil@pmeasuring.com

CHINA
T: +86 400 081 8020
E: pmschina@pmeasuring.com

FRANCE
T: +33 682 991 798
E: pmsfrance@pmeasuring.com

GERMANY
T: +49 351 8896 3850
E: pmsgermany@pmeasuring.com

IRELAND
T: +353 1 295 7373
E: Info.ie@pmeasuring.com

ITALY
T: +39 06 9053 0130
E: pmsrl@pmeasuring.com

JAPAN
T: +81 44 589 3498
E: pmsjapan@pmeasuring.com

KOREA
T: +82 31 286 5790
E: pmskorea@pmeasuring.com

NORDIC
T: +45 7070 2855
E: pmsnordic@pmeasuring.com

PUERTO RICO
T: +1 787 718 9096
E: pmspuertorico@pmeasuring.com

SINGAPORE
T: +65 6496 0342
E: pmsingapore@pmeasuring.com

SWITZERLAND
T: +41 71 987 01 01
E: pmsswitzerland@pmeasuring.com

TAIWAN
T: 886-3-5525300 Ext: 301
E: pms taiwan@pmeasuring.com

UNITED KINGDOM
T: +44 (0)1223 257704
E: Info.uk@pmeasuring.com

For the most updated contacted information please visit <https://www.pmeasuring.com/contact-us/>



LiQuilaz®, Liquistat®, NanoVision Technology®, UltraChem®, and Ultra DI® are registered trademarks of Particle Measuring Systems®, Inc. All other trademarks are the property of their respective owners. Particle Measuring Systems, Inc. reserves the right to change specifications without notice. © 2024 Particle Measuring Systems, Inc. All rights reserved. 09172024